


New opportunities for paramagnetic materials using ultra-fast MAS

Frank Engelke, Sebastian Wegner

Bruker BioSpin GmbH

EU-FP7: pNMR Workshop, Aarhus, 8 July 2015

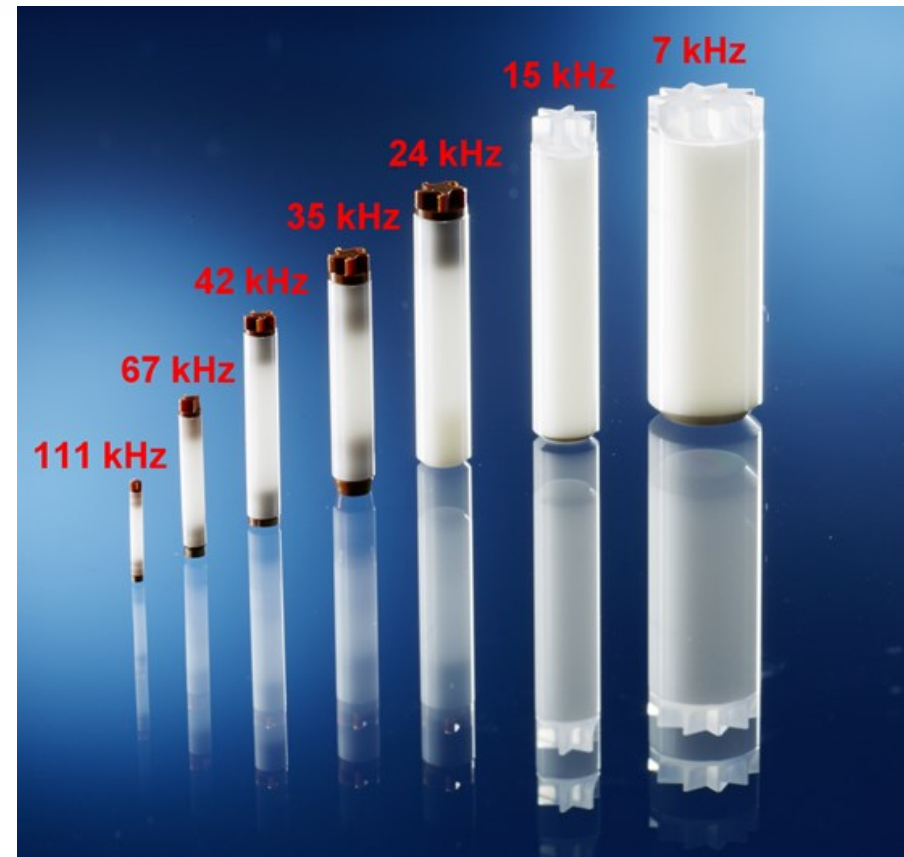

Collaborations and acknowledgements



BBIO

Benno Knott
David Osen
David Bennett (ER)
Sebastian Wegner
Armin Porea
Andreas Ruppel
Alexander Krahm
Heiko Beiler
Reinhard Steppe
Fabien Aussenac

ENS Lyon

Guido Pintacuda
Loren Andreas
Daniela Lalli
Kevin Sanders
Anne Lesage
Sachin Chaudhuri

ZHAW Winterthur

Dirk Wilhelm
Nicoleta Herzog
Alexander S. Weber



Outline

1. Why faster MAS ?
2. What means fast MAS ?
3. Probe hardware for fast MAS
4. How about CRAMPS ?
5. Heteronuclear decoupling under fast MAS
6. Cross polarization under fast MAS
7. Some applications to paramagnetic systems
8. Fast MAS and extended VT

Why faster and faster Rotation ?

- Strong dipole – dipole interactions
 - Homonuclear dipole – dipole couplings
 - $^1\text{H} \leftrightarrow ^1\text{H}$ and $^{19}\text{F} \leftrightarrow ^{19}\text{F}$ (also $^1\text{H} \leftrightarrow ^{19}\text{F}$)
 - Heteronuclear dipole – dipole couplings
 - MAS frequency should be 3-5 times higher than the coupling
- Quadrupolar nuclei
 - Fast MAS rotation enables the observation of undistorted central transitions, free of spinning sidebands
 - Spinning sidebands from satellites can be detected more easily undistorted
- Paramagnetic species
 - Paramagnetically distorted samples can have anisotropies resulting in spectra with widths more than a 1.000 ppm

Some common interactions in solid state NMR




Parameter	Order of magnitude	Example
CSA, chemical shift	up to 100 kHz	^{77}Se , ^{119}Hg
Dipole dipole coupling	up to 100 kHz	^{19}F , (^1H)
J-coupling	up to 10 kHz	$^1\text{J}(^{31}\text{P}, ^{31}\text{P})$
Quadrupolar coupling	If eQ small, only 1 st order important, can be averaged by MAS	^2H , ^{14}N , ^6Li
	If eQ large, than 1 st and 2 nd order are important, second order not averaged but only scaled by MAS	^{27}Al $^{47/49}\text{Ti}$

-> fast MAS in the regime of 100kHz can be beneficial for some interactions !

What means fast MAS?

The terms “fast MAS” or “very fast MAS” or “ultrafast MAS”
are a function $f(t)$ of time t .

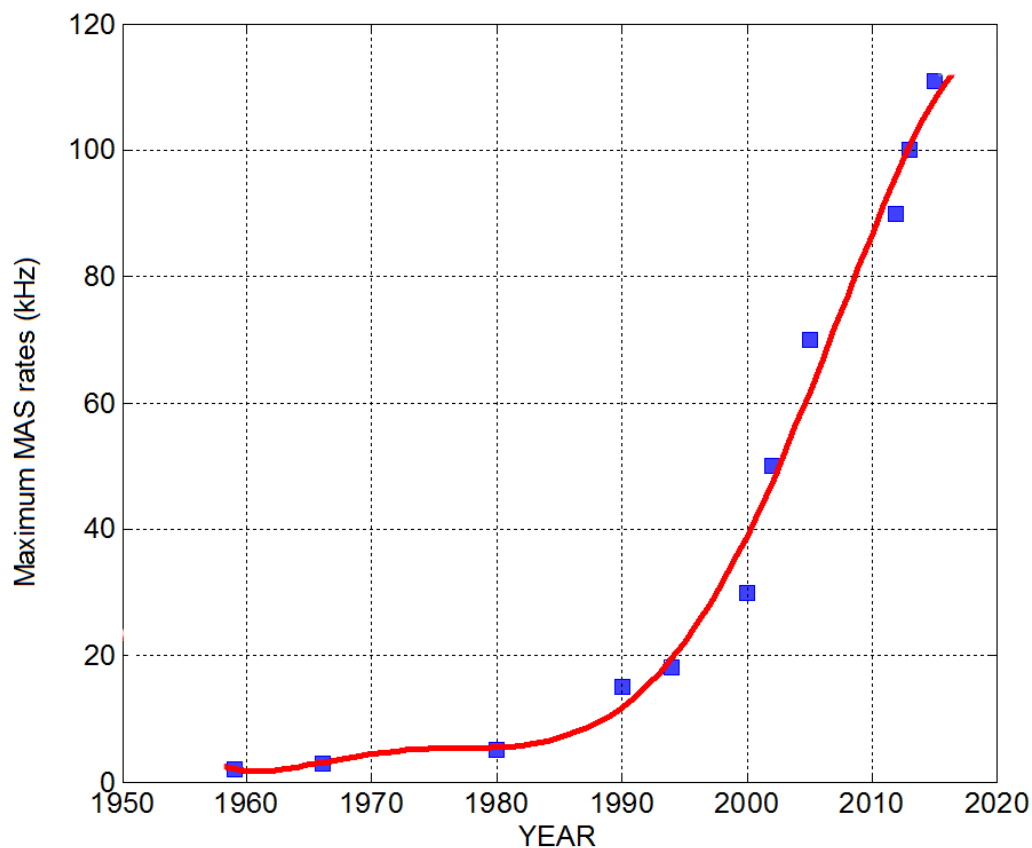

Technology for fast MAS


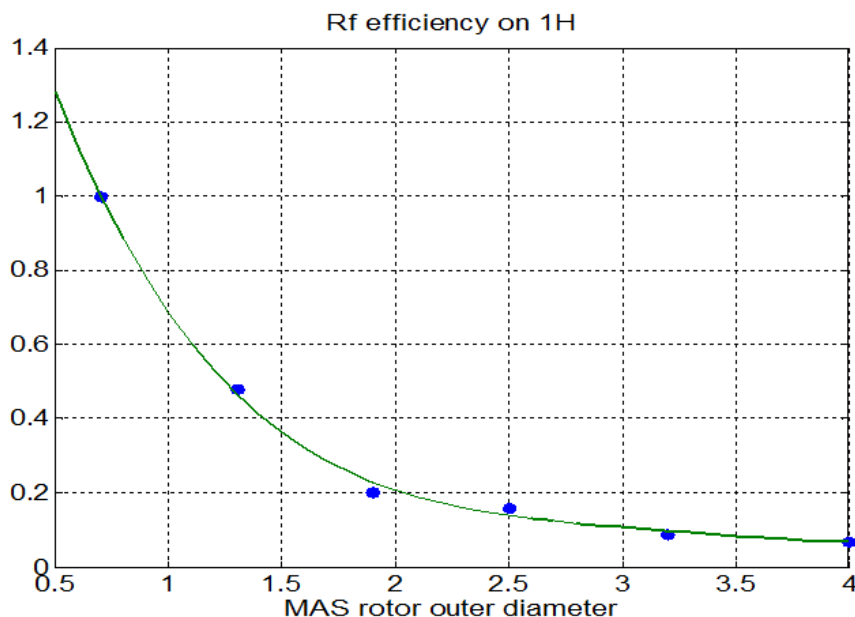


Fig. 7: Relative ^1H rf efficiencies for MAS probes with rotor diameters from 0.7mm to 4mm.

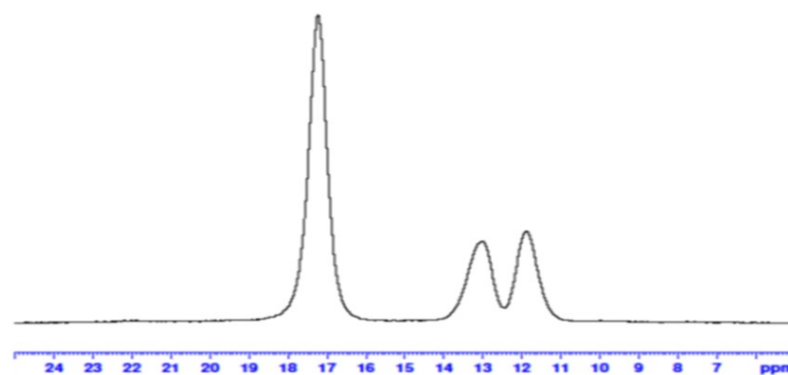

^1H spectrum of glycine at 1000 MHz and spinning at an MAS speed of 111 kHz.

Single pulse excitation
($P1 = 1.1\mu\text{s}$), single scan.

Active sample volume 350 nl

MAS 0.7mm


Control of bearing and drive flow

Frictional heating



MAS	CS difference	frictional heating	sample T
10 kHz	0	0	15
40 kHz	0	0	15
60 kHz	0,1	4	19
80 kHz	0,3	12	27
100 kHz	0,5	20	35
111 kHz	0,6	24	39
111 kHz + BCU-X	0	0	15


**MAS 3 PNEUMATIC
CONTROL UNIT**

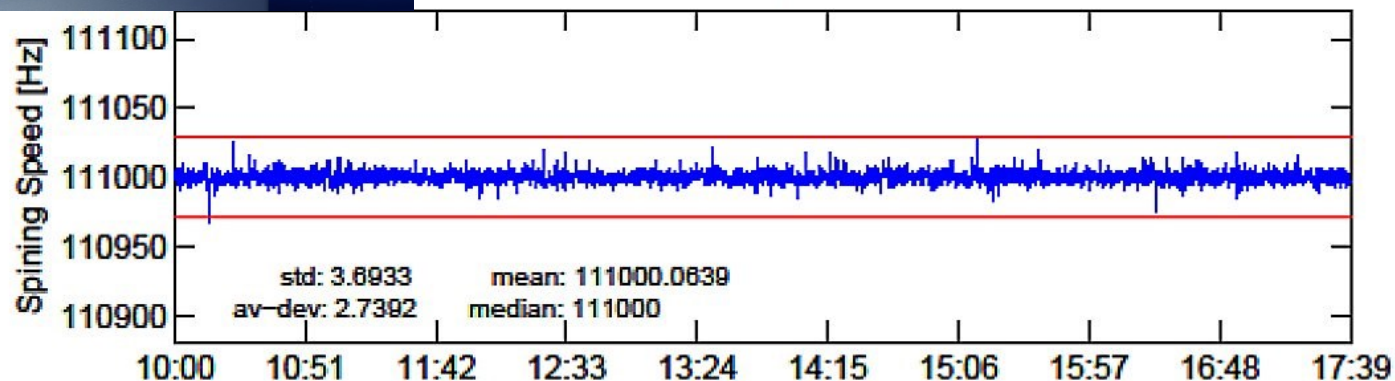


Technology for fast MAS – MAS 3

- Spin regulation up to 200kHz
- Fiberoptical spin-signal-interface to the probe
(and backward compatible spin-interface)
- New spinrate readout resolution: 0,1Hz
- Additional pneumatic channels for future probe designs
- Safer and faster communication using ethernet and CAN-bus
- Several logging functions
- Internal 3-axis magnetic field sensors for coarse stray field measurements
- Easy probe profile management therefore easy A2B conversion

Fast MAS – technological limits

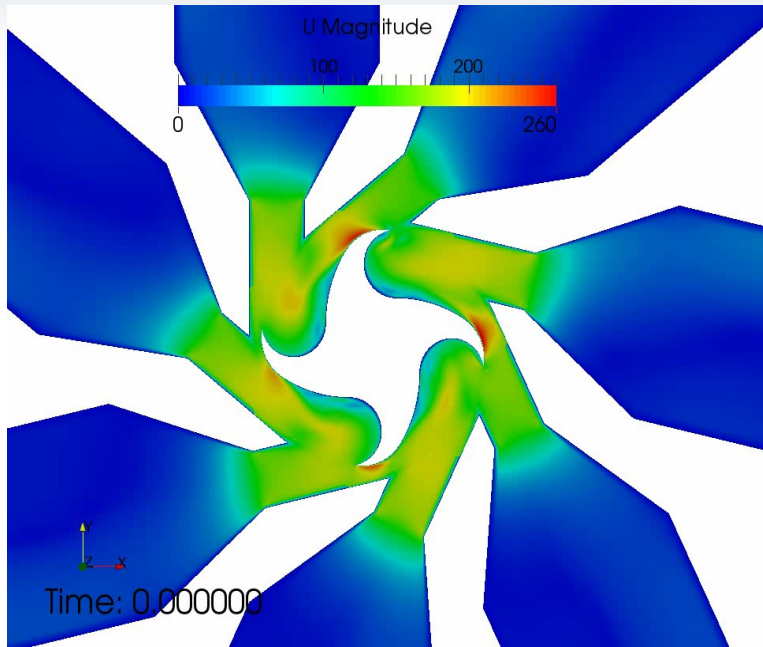


Gas velocity profiles for
MAS 1.3mm turbines

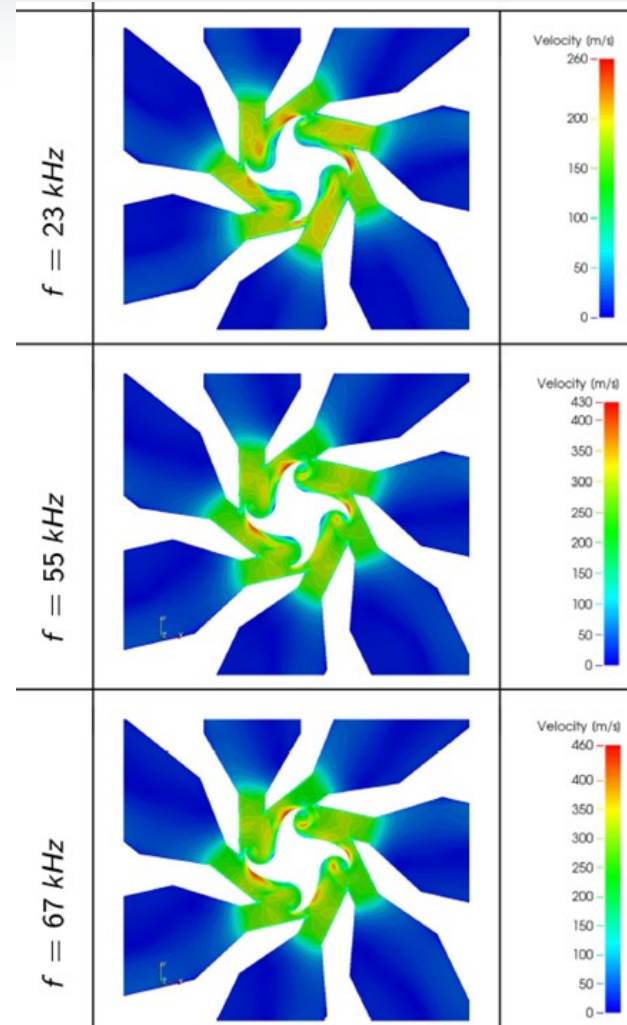

Spectral resolution: What about the good old CRAMPS ?

- @ 70kHz -> 1.3mm probe
the result seems
disappointing, since
70kHz is not fast
enough

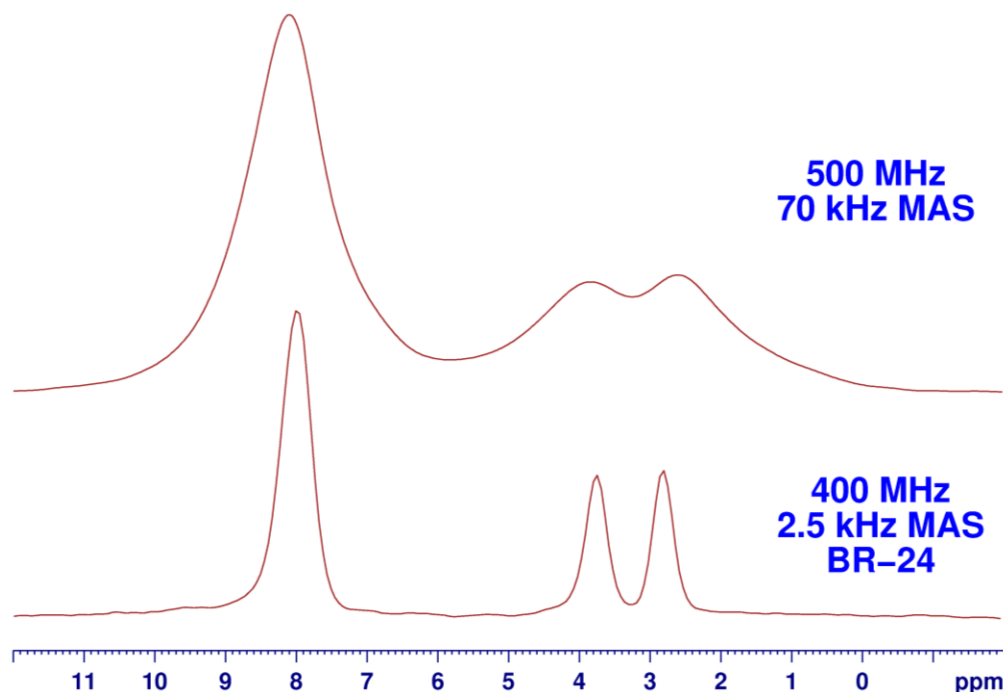


Resolution: What about the good old CRAMPS ?

- @ 70kHz → 1.3mm probe the result seems to be insufficient, since spinning at 70kHz is not fast enough.

- @ 111kHz it looks better

→ 111kHz exceeds the homonuclear coupling strength of glycine

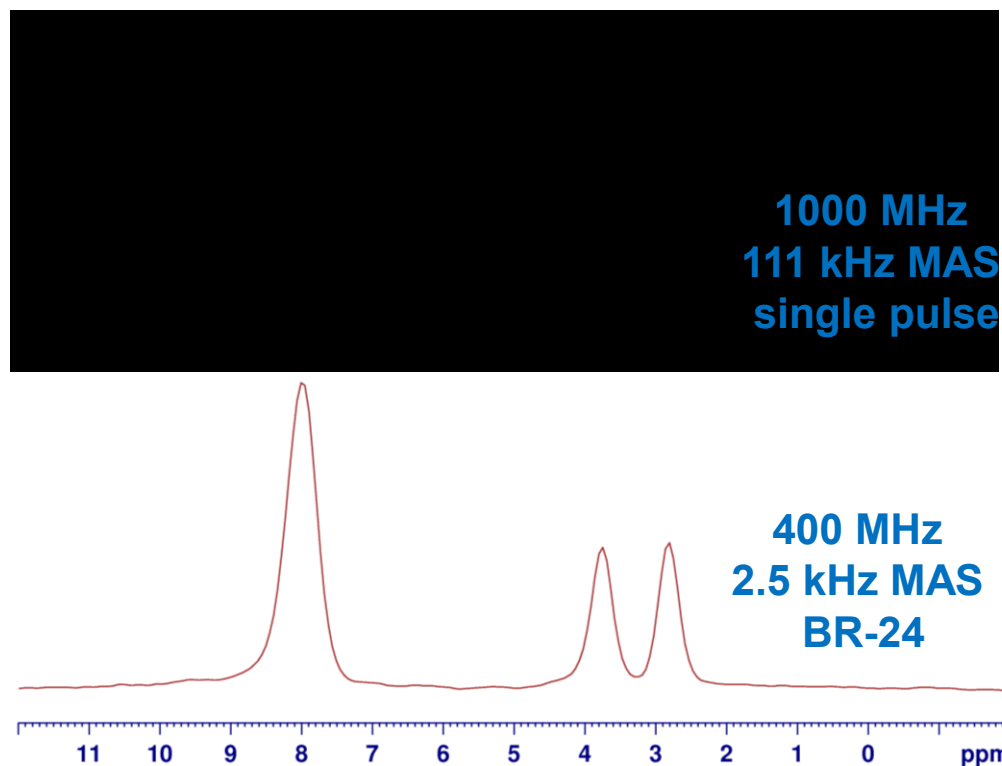


Decoupling under fast MAS




CW Decoupling

- Heteronuclear decoupling can be done with different approaches
 - High power decoupling
 - Low power CW decoupling matched to MAS speed

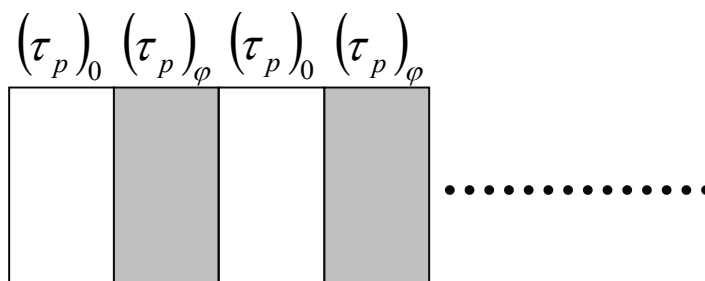

TPPM / SPINAL

CW Decoupling at 42kHz (MAS 1.9mm)

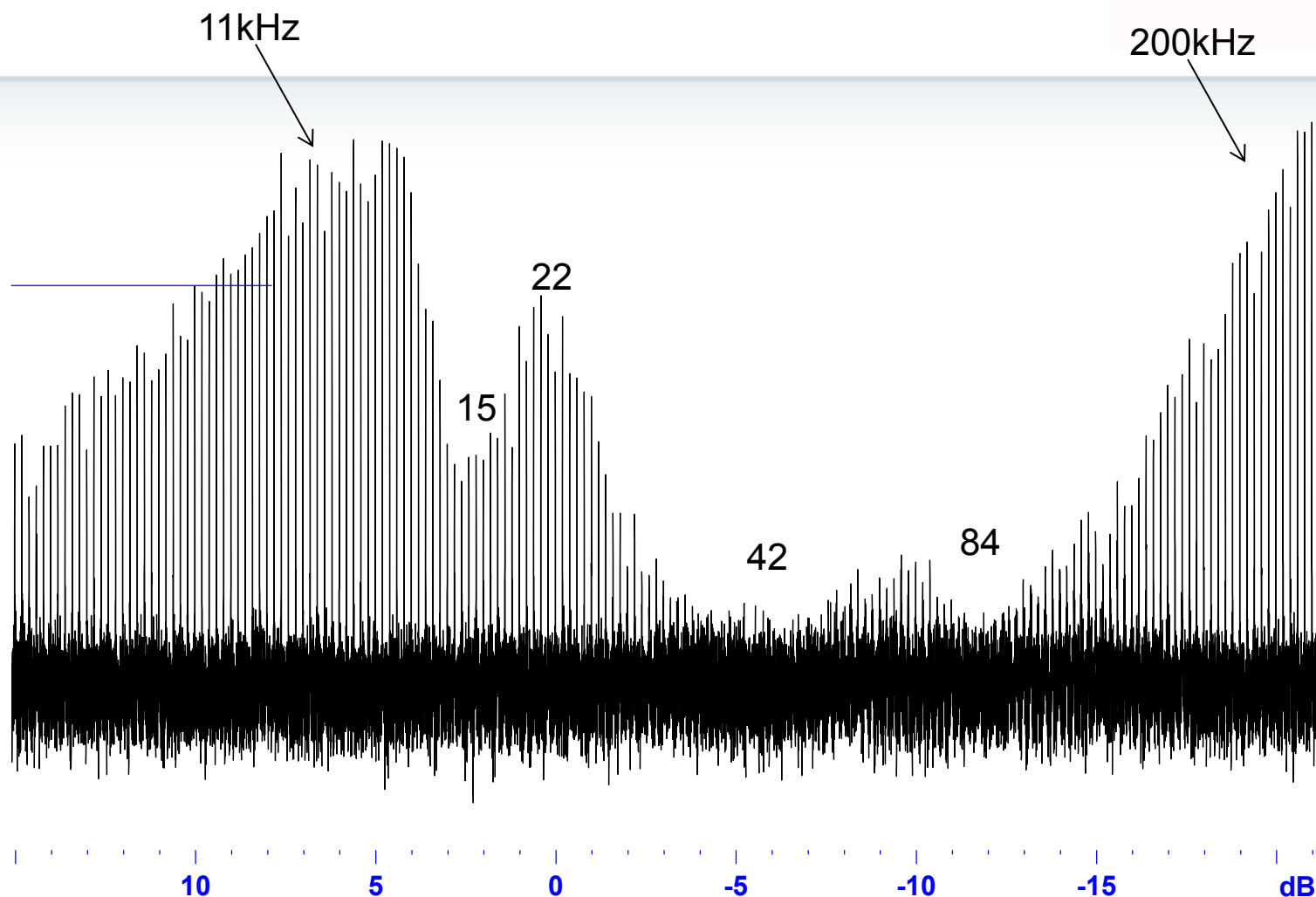


See, e.g., M. Weingarth, et al., Chem.Phys.Lett. 466, 247-251, 2009.
M. Ernst, J. Magn. Reson. 162, 1-34, 2003.

CP efficiency at fast MAS

example @ 42kHz (MAS 1.9mm)




- Approaching very fast Magic Angle Spinning the dipolar coupling gets averaged -> what about cp efficiencies?
- Intensity of $^{13}\text{CH}_2$ signal in glycine at various spinning rates
 - Same contact time
 - Same ^{13}C RF-field
 - ^1H RF-field was adjusted while (using a 50% to 100% ramp)
 - MAS 10 kHz
 - MAS 25 kHz
 - MAS 42 kHz

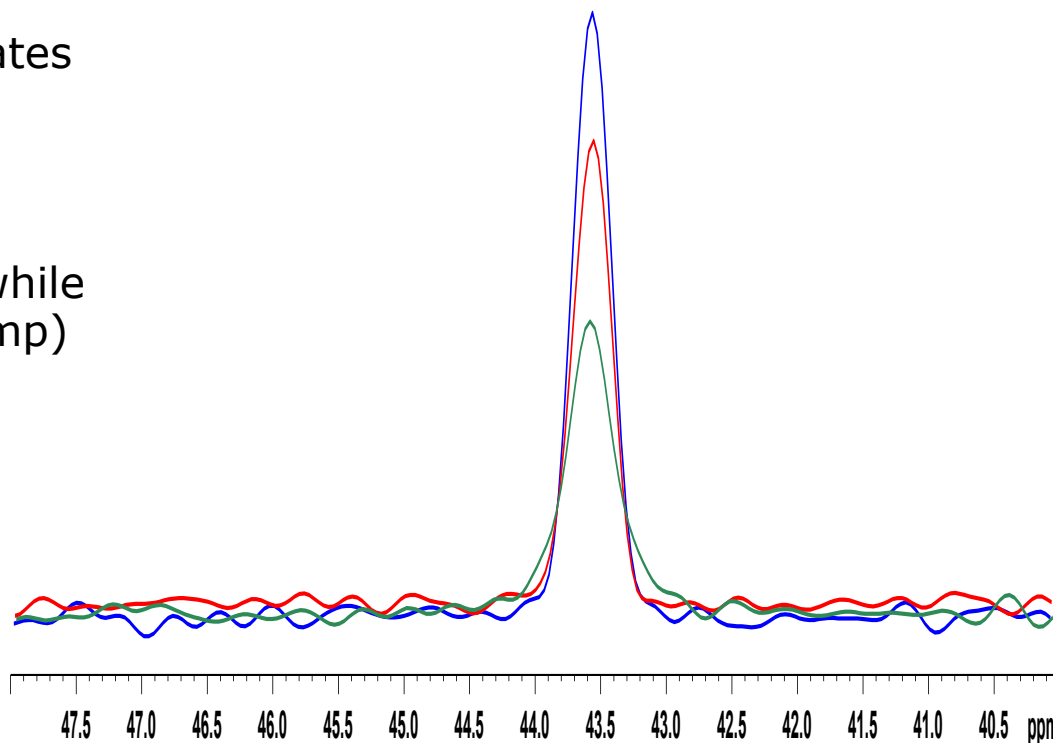


CP and MAS




(a) Broadening of energy levels due to I - I dipolar couplings

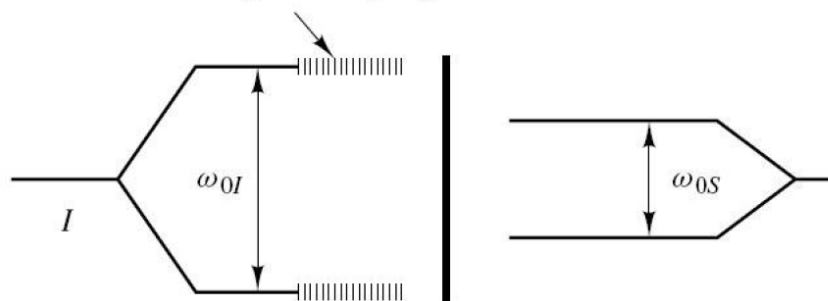

(b)

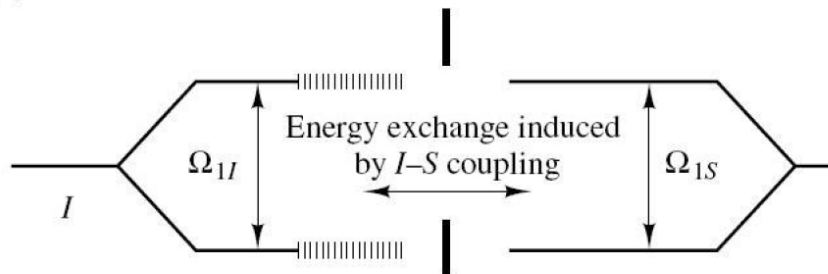

Hartmann-Hahn matching: $\Omega_{1I} = \Omega_{1S}$

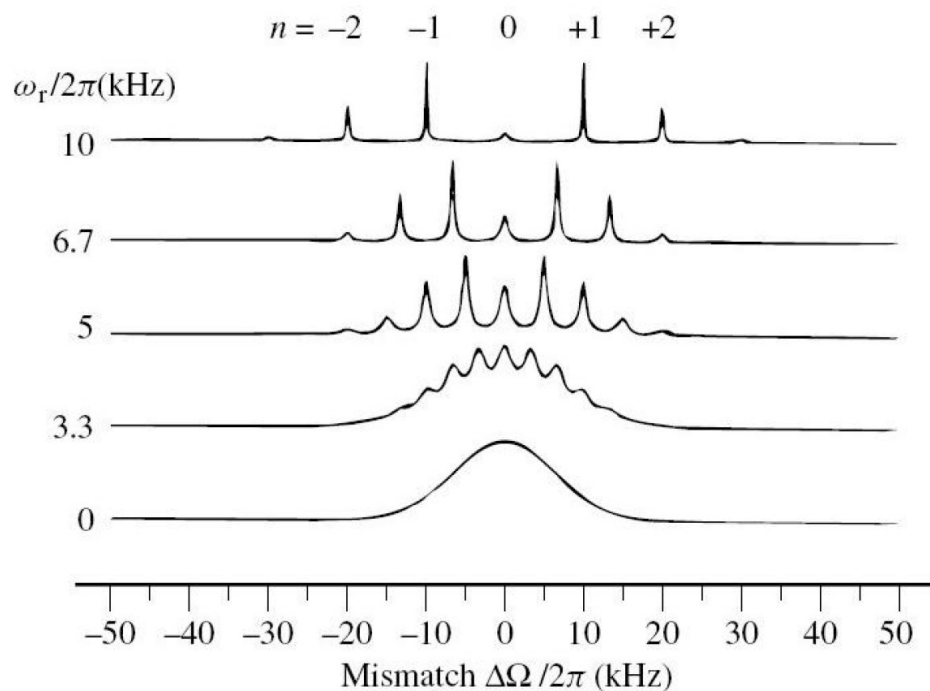

See, e.g., E.O. Stejskal, et al, J. Magn. Reson. 57, 471, 1984

CP MAS dynamics




Using MAS the energy levels in the rotating frame are modulated by the magic angle rotation.

For the case when ω_R is small compared to the RF-fields $\omega_{1,I}$ and $\omega_{1,S}$ the CP Hartmann-Hahn matching conditions is given by:

$$\omega_{1,I} \pm n \omega_R = \omega_{1,S}$$

$$\omega_{1,I} - \omega_{1,S} = \pm n \omega_R$$

This is shown on the right side for $n=1$: The polarisation transfer is supported by *flip-flop* terms (zero-quantum transitions)

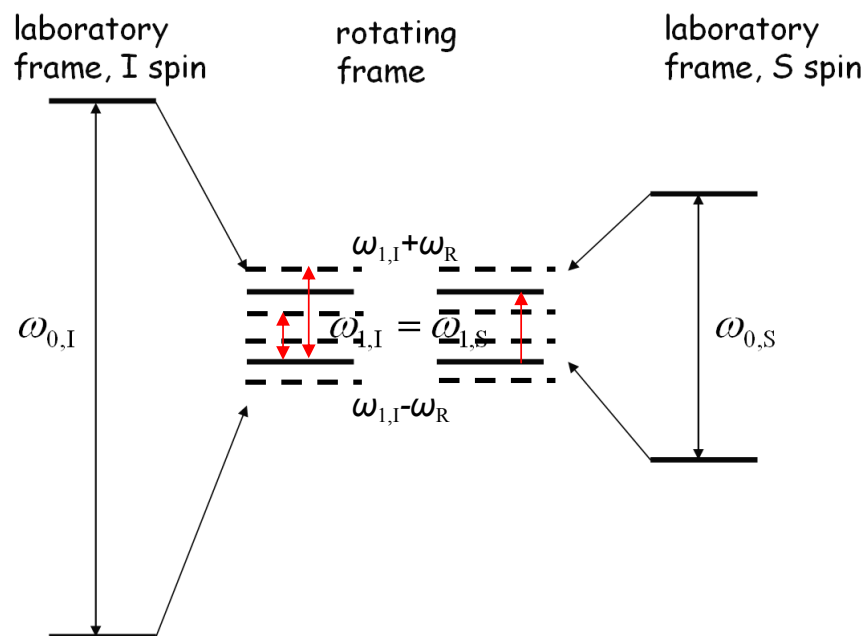

CP and very fast MAS




When the MAS speed ω_R is approaching a regime, where ω_R is faster (stronger) than the RF-fields $\omega_{1,I}$ and $\omega_{1,S}$ the CP Hartmann-Hahn condition can also be:

$$\omega_{1,I} \pm n \omega_R = -\omega_{1,S}$$

$$\omega_{1,I} + \omega_{1,S} = \pm n \omega_R$$

Shown on the right for $n = 1$; now the polarisation transfer is also supported by *flop-flop* processes, the elemental CP step is now a double-quantum transition.

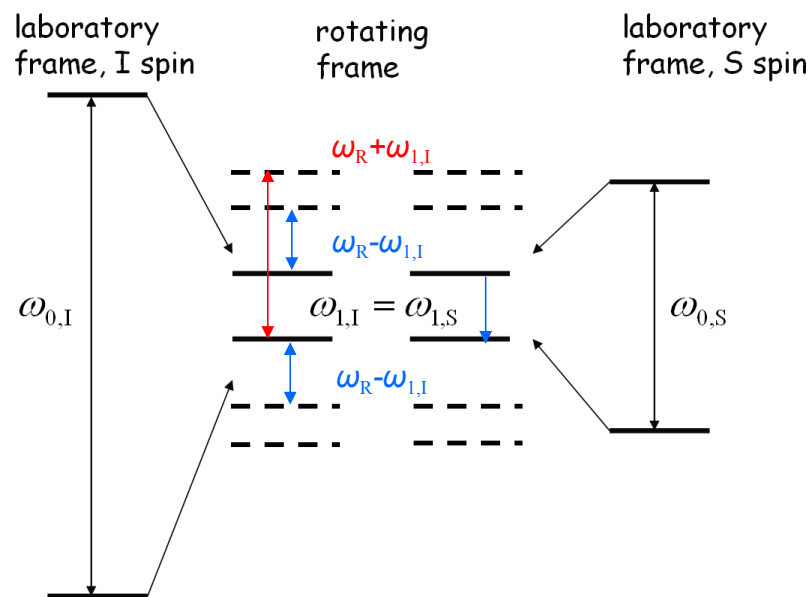

See, e.g., S. Laage, et al., J. Magn. Reson. 196, 133, 2009

Hartmann-Hahn at 111kHz MAS

ZQ condition:

$$\frac{\omega_{1I} - \omega_{1S}}{\omega_R} = \pm n$$


DQ condition:

$$\frac{\omega_{1I} + \omega_{1S}}{\omega_R} = \pm n$$

For $\omega_R = 111$ kHz, $\omega_{1S} = 15$ kHz
matching conditions:

$\omega_{1I}(\text{ZQ}, n=0) = 15$ kHz
 $\omega_{1I}(\text{ZQ}, n=1) = 126$ kHz
 $\omega_{1I}(\text{ZQ}, n=2) = 237$ kHz
 $\omega_{1I}(\text{ZQ}, n=3) = 348$ kHz

$\omega_{1I}(\text{DQ}, n=1) = 96$ kHz
 $\omega_{1I}(\text{DQ}, n=2) = 207$ kHz
 $\omega_{1I}(\text{DQ}, n=3) = 318$ kHz

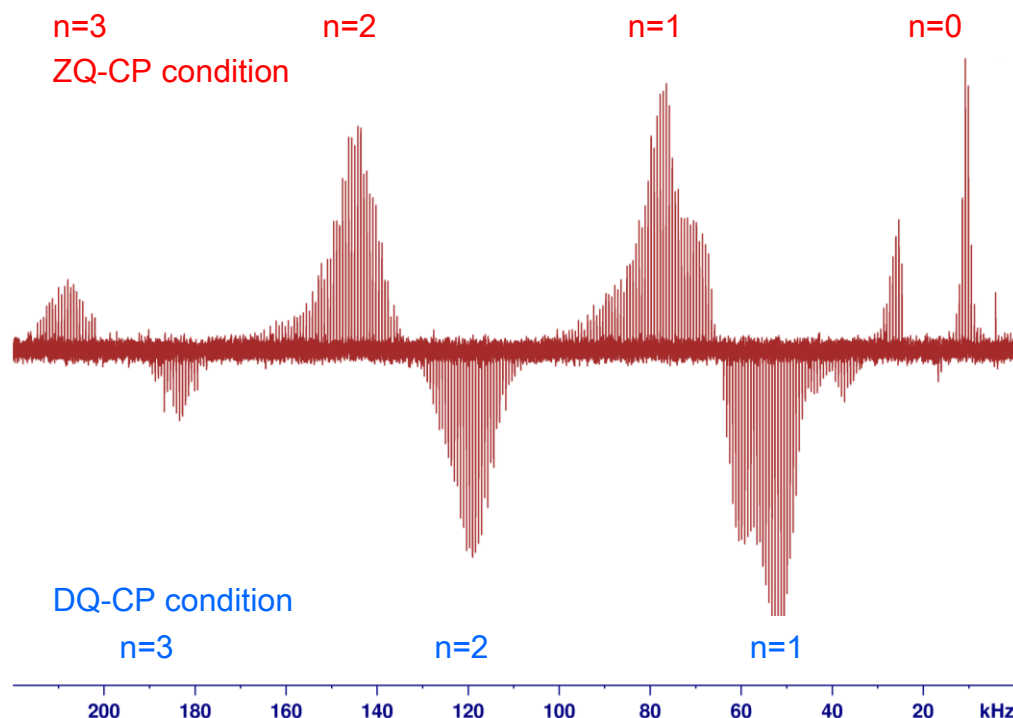

Hartmann-Hahn at 111kHz MAS destructive overlay of H-H conditions

ZQ Kondition:

$$\frac{\omega_{1S} - \omega_{2S}}{\omega_R} = \pm n$$


DQ Kondition:

$$\frac{\omega_{1S} + \omega_{2S}}{\omega_R} = \pm n$$

Für $\omega_R = 111$ kHz, $\omega_{1I} = 55$ kHz
the matching conditions are:

$$\left. \begin{aligned} \omega_{1S}(\text{ZQ}, n=0) &= 55 \text{ kHz} \\ \omega_{1S}(\text{ZQ}, n=1) &= 166 \text{ kHz} \\ \omega_{1S}(\text{ZQ}, n=2) &= 277 \text{ kHz} \\ \omega_{1S}(\text{ZQ}, n=3) &= 388 \text{ kHz} \end{aligned} \right\}$$

$$\left. \begin{aligned} \omega_{1S}(\text{DQ}, n=1) &= 56 \text{ kHz} \\ \omega_{1S}(\text{DQ}, n=2) &= 167 \text{ kHz} \\ \omega_{1S}(\text{DQ}, n=3) &= 278 \text{ kHz} \end{aligned} \right\}$$


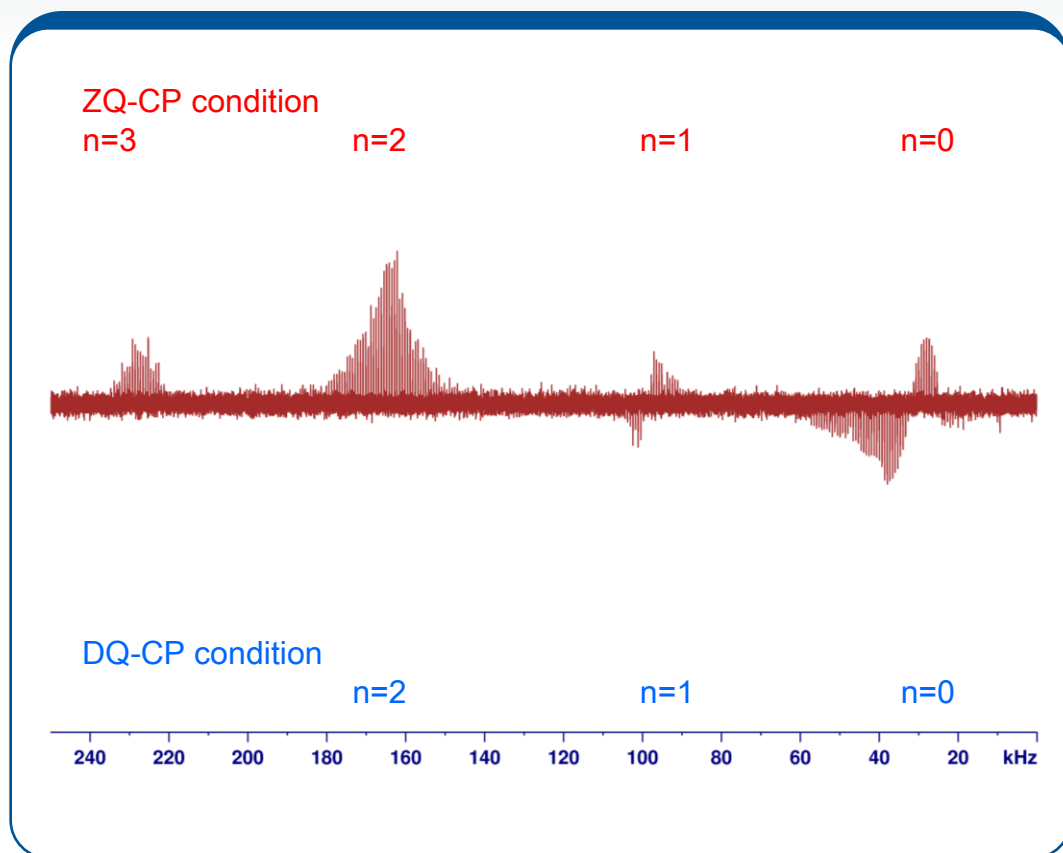


MAS 0.7 mm

^1H detected ^{15}N - ^1H 2D correlation @ 111kHz




Coherence Lifetimes & Linewidths (GB1)

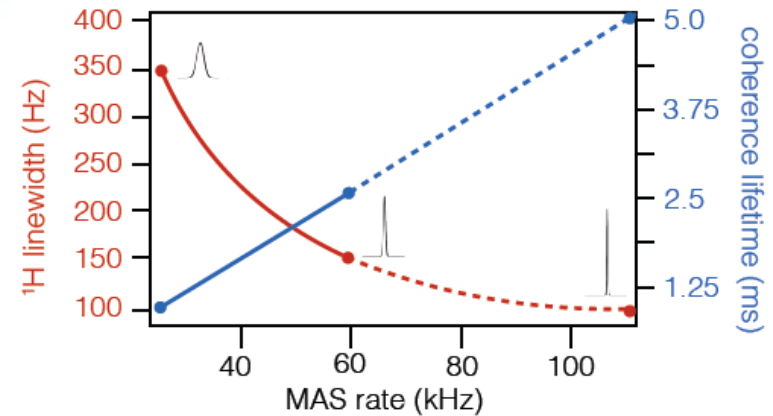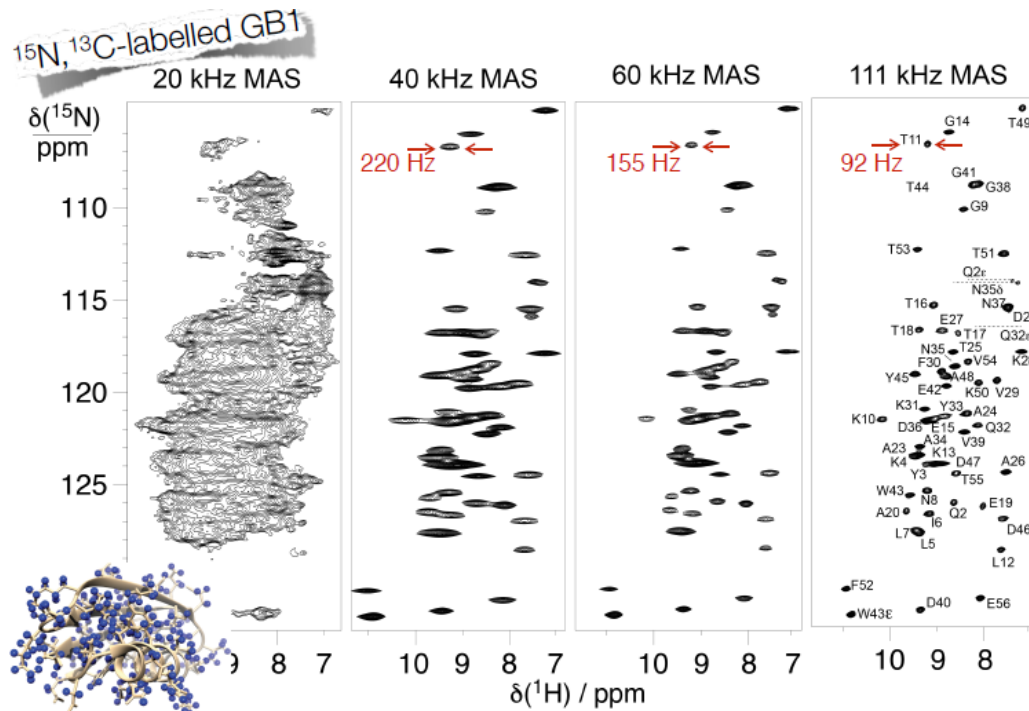



Paramagnetic metalloproteins: Metal-ion coordination



$(^{13}\text{C}, ^{15}\text{N})$ HiPIP 1 from *E. halophila*




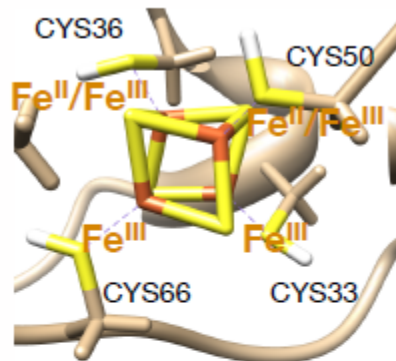


$(^2\text{H}, ^{13}\text{C}, ^{15}\text{N})$ $\text{Cu}^{\text{II}}, \text{Co}^{\text{II}}$ -human superoxide dismutase


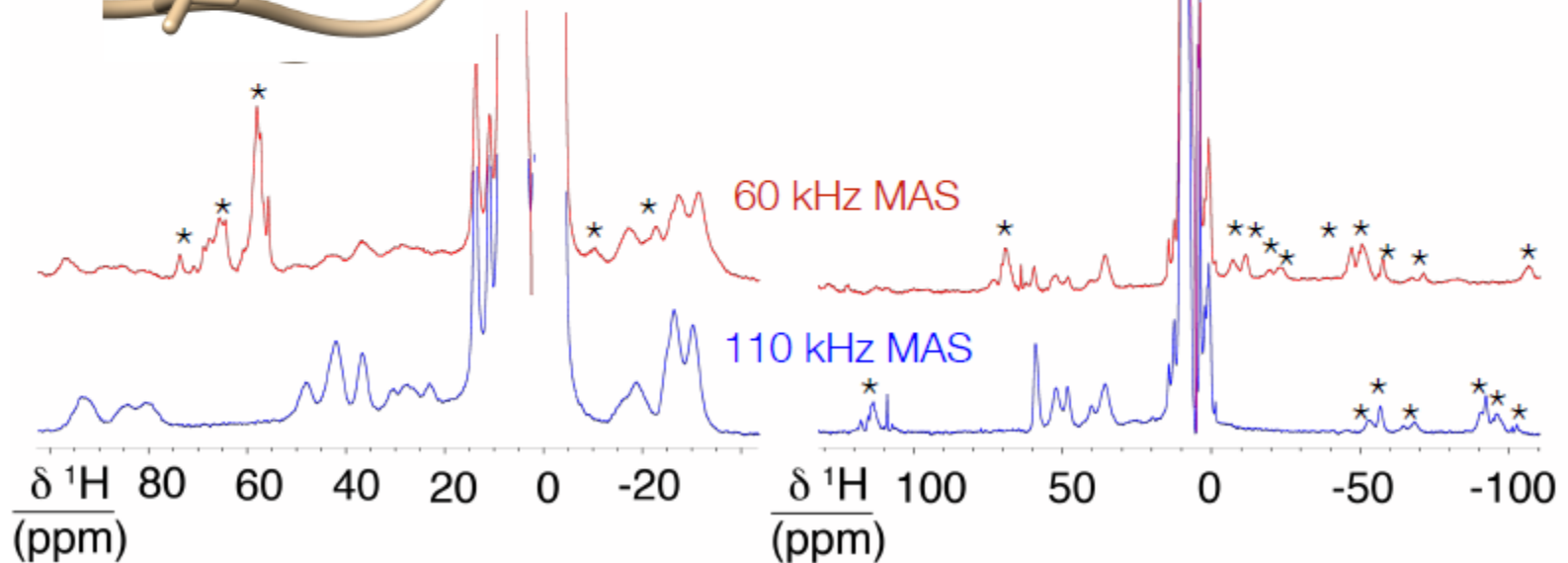
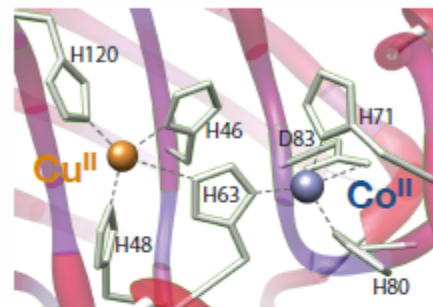


Paramagnetic Solids

Large Offsets and Large Shift Anisotropies

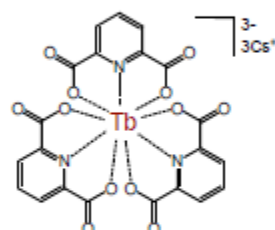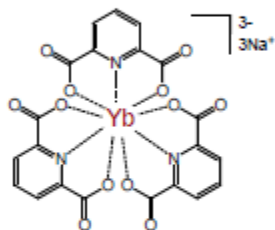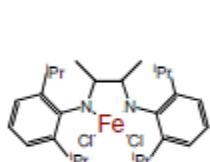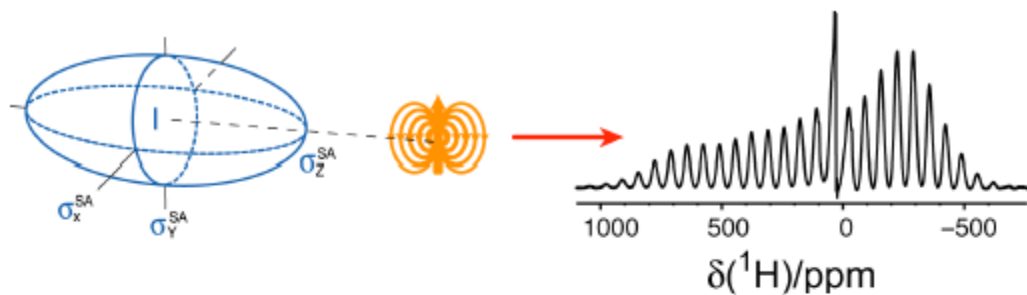

pNMR Network

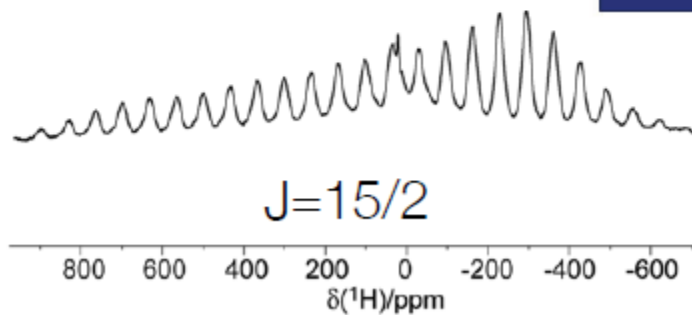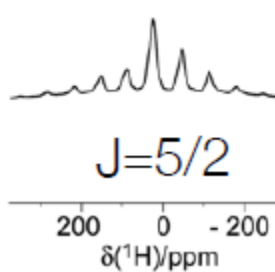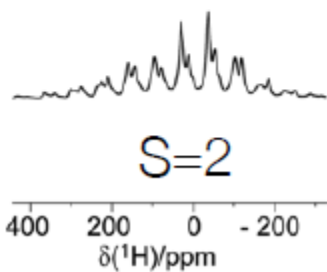

Fast MAS and extended VT

- Ultrafast MAS (0.7 mm)
(ENS Lyon February 2015)
- Fast MAS (1.3 mm) low temperature (100 K) DNP probe
(ENS Lyon 2016, ETHZ 2017)
- Fast MAS (1.3 mm) non-DNP with extended VT capabilities
(functional prototype July 2016)

*Thank you
for your attention !*